

Cubis[®] MCM5004

Manual Mass Comparator

User Benefits

- Complete mass standard laboratory in a single unit
- Integrated climate sensors for recording all data relevant for determining measurement uncertainty
- Integrated workflow control for efficient and error-free mass comparison
- Fast measurement cycles according to the ABA, ABBA or AB₁...B_nA method



Highlighted Performance Features

- Cubis[®] MSA color touch screen for fast and simple configuration of parameters and workflows
- Sensor-equipped climate module integrated into the draft shield for recording the temperature, humidity and air pressure
- Integrated calibration workflows for ABA, ABBA, AB₁...B_nA cycles to ensure efficient, error-free mass comparison
- Fully integrated function for determining the measurement uncertainty in accordance with OIML and ASTM recommendations
- Filters for optimal adaptation of the mass comparator to ambient conditions
- Monolithic weighing technology
- For display and evaluation, complete electronics and power supply separated from the weighing system to prevent heat from affecting the results
- All MCM mass comparators featuring eccentric (off-center) load compensation for easy loading of weights without automatic centering
- Additional applications for density determination, statistics and individual identifiers are integrated as standard programs
- Automatic, motorized leveling
- Built-in SD card slot for storage and transfer of all data and settings
- Easy logging of reference weight data
- Continuous weighing range display: any weight between 0 g and the maximum capacity can be displayed
- USB, RS-232C and Ethernet interface ports to integrate the mass comparator into networks or to enable it to communicate with external software via third-party protocols, standardized communication protocols or Web services

Technical Specifications

Metrological Specifications	
Maximum capacity	5100 g
Application range	0 – 5100 g
Readability	0.1 mg
Repeatability, optimal ¹⁾	0.3 mg
Repeatability, standard E ²⁾	0.5 mg
Repeatability, E ¹ / ₁₀ load ²⁾	0.3 mg
Repeatability standard, F ³⁾	0.8 mg
Electronic weighing range and tare range	5100 g
Linearity	2 mg
Eccentric load deviation	151 µg mm
Stabilization time	3 s
Cycle time, ABBA in s	90 s

Basic Equipment	
Interfaces	RS232C USB LAN
isoCAL	✓
Draft shield	✓
Application programs	Basic weighing, mass unit conversion, individual identifiers, density determination, statistics
Below-comparator weighing port	✓
Air temperature sensor	✓
Air humidity sensor	✓
Air pressure sensor	✓
PC connecting cable	USB

Ambient Conditions	
Permissible operating temperature range	10–30 °C
Recommended operating temperature	22 °C
Temperature fluctuations	0.3°C/h 0.5°C/12h
Max. air current	< 0.2 m/s
Humidity range	40 – 70 %
Humidity fluctuations	5% 4 h
Power supply	100 – 240 V AC/50 – 60 Hz
Power consumption	< 35 VA

Dimensions	
Weighing pan diameter	136 × 136 mm
Sample size (D × H)	130 × 200 mm
Weigh cell (W × D × H)	240 × 276 × 373 mm
Electronic unit (W × D × H)	239 × 320 × 56 mm
Net weight	15 kg
Gross weight	22.5 kg
Number of packages	1
Packaging data 1	83 × 45 × 59 cm
Optimal height for setup	800 mm

Applications	
OIML R111, class E1	5 kg
OIML R111, class E2	2 – 5 kg
OIML R111, class F1	200 g – 5 kg
OIML R111, class F2	100 g – 5 kg
OIML R111, class M1	
OIML R111, class M2	
OIML R111, class M3	
ASTM E617, class 0	1 – 5 kg
ASTM E617, class 1	500 g – 5 kg
ASTM E617, class 2	300 g – 5 kg
ASTM E617, class 3	100 g – 5 kg
ASTM E617, class 4	
ASTM E617, class 5	
ASTM E617, class 6	

Optional Accessories	
External calibration weight	5 kg E2 YCW652-00
Climate module, uncalibrated, for all MCM models	YCM20MC
Calibration of a YCM20MC climate module with DAkkS calibration certificate	YCM20DAkkS
Climate module with DAkkS calibration certificate for all MCM models	YCM20MC-DAkkS
Optional draft shield	YDS24C
Weighing table	YWT03

The standard deviation "s" is the repeatability calculated from 5 ABA cycles under the following conditions:

- ¹⁾ Optimal conditions: automatic measurement without operator influence measured in a laboratory under E1 conditions, on a decoupled weighing stone no drafts from above
- ²⁾ Standard conditions E: measured by hand in a laboratory under E1 conditions, on a decoupled weighing stone; no drafts from above
- ³⁾ Standard conditions F: measurement performed manually in a laboratory under at least F1 conditions, on a non-decoupled weighing stone, air conditioning and minimal drafts from above

Sales and Service Contacts

For further contacts, visit www.sartorius.com

Europe

Germany

Sartorius Lab Instruments
GmbH & Co. KG
Weender Landstrasse 94-108
37075 Goettingen

Phone +49.551.308.0
Fax +49.551.308.3289

France & Suisse Romande

Sartorius France
2, rue Antoine Laurent de Lavoisier
ZA de la Gaudrée
91410 Dourdan

Phone +33.1.70.62.50.00
Fax +33.1.64.59.76.39

Austria

Sartorius Austria GmbH
Franzosengraben 12
1030 Vienna

Phone +43.1.7965760.0
Fax +43.1.7965760.24

Belgium

Sartorius Belgium N.V.
Leuvensesteenweg, 248/B
1800 Vilvoorde

Phone +32.2.756.06.71
Fax +32.2.253.45.95

Finland & Baltics

Sartorius Biohit Liquid Handling Oy
Laippatie 1
00880 Helsinki

Phone +358.9.755.951
Fax +358.9.755.95.200

Hungary

Sartorius Hungária Kft.
Kagyló u. 5.
2092 Budakeszi

Phone +3623.457.227
Fax +3623.457.147

Ireland

Sartorius Ireland Ltd.
Unit 41, The Business Centre
Stadium Business Park
Ballycoolin Road
Dublin 11

Phone +353.1.8089050
Fax +353.1.8089388

Italy

Sartorius Italy S.r.l.
Viale A. Casati, 4
20835 Muggiò (MB)

Phone +39.039.4659.1
Fax +39.039.4659.88

Netherlands

Sartorius Netherlands B.V.

Phone +31.30.60.53.001
Fax +31.30.60.52.917
info.netherlands@sartorius.com

Poland

Sartorius Poland sp. z o.o.
ul. Wrzesinska 70
62-025 Kostrzyn

Phone +48.61.6473830
Fax +48.61.6473839

Russian Federation & Ukraine

LLC "Sartorius ICR"
and LLC "Biohit"
Uralskaya str. 4, Lit. B
199155, Saint-Petersburg

Phone +7.812.327.5.327
Fax +7.812.327.5.323

LLC "Biohit"
Post Box 440 "B"
01001 Kiev, Ukraine

Phone +380.44.411.4918
Fax +380.50.623.3162

Spain & Portugal

Sartorius Spain, S.A.
Avda. de la Industria, 32
Edificio PAYMA
28108 Alcobendas (Madrid)

Phone Spain +34.902.123.367
Phone Portugal +351.800.855.800
Fax Spain +34.91.358.96.23
Fax Portugal +351.800.855.799

Switzerland

Sartorius Mechatronics Switzerland AG
Ringstrasse 24a
8317 Tagelswangen (ZH)

Phone +41.44.746.50.00
Fax +41.44.746.50.50

U.K.

Sartorius UK Ltd.
Longmead Business Centre
Blenheim Road, Epsom
Surrey KT19 9QQ

Phone +44.1372.737159
Fax +44.1372.726171

America

USA

Sartorius Corporation
5 Orville Drive, Suite 200
Bohemia, NY 11716

Phone +1.631.254.4249
Toll-free +1.800.635.2906
Fax +1.631.254.4253

Argentina

Sartorius Argentina S.A.
Int. A. Ávalos 4251
B1605ECS Munro
Buenos Aires

Phone +54.11.4721.0505
Fax +54.11.4762.2333

Brazil

Sartorius do Brasil Ltda
Avenida Senador Vergueiro 2962
São Bernardo do Campo
CEP 09600-000 - SP- Brasil

Phone +55.11.4362.8900
Fax + 55.11.4362.8901

Canada

Sartorius Canada Inc.
2179 Dunwin Drive #4
Mississauga, ON L5L 1X2

Phone +1.905.569.7977
Toll-Free +1.800.668.4234
Fax +1.905.569.7021

Mexico

Sartorius de México S.A. de C.V.
Circuito Circunvalación Poniente
No. 149
Ciudad Satélite
53100, Estado de México
México

Phone +52.5555.62.1102
Fax +52.5555.62.2942

Asia | Pacific

Australia

Sartorius Australia Pty. Ltd.
Unit 5, 7-11 Rodeo Drive
Dandenong South Vic 3175

Phone +61.3.8762.1800
Fax +61.3.8762.1828

China

Sartorius Scientific
Instruments (Beijing) Co., Ltd.
33 Yu An Road, Airport Industrial Park
Zone B,
Shunyi District, Beijing 101300,
P.R.China

Phone +86.10.8042.6300
Fax +86.10.8042.6486

Hong Kong

Sartorius Hong Kong Ltd.
Unit 1012, Lu Plaza
2 Wing Yip Street
Kwung Tong
Kowloon, Hong Kong

Phone +852.2774.2678
Fax +852.2766.3526

India

Sartorius Weighing India Pvt. Ltd.
#69/2-69/3, NH 48, Jakkasandra,
Nelamangala Tq
562 123 Bangalore, India

Phone +91.80.4350.5250
Fax +91.80.4350.5253

Japan

Sartorius Stedim Japan K.K.
4th Fl., Daiwa Shinagawa North Bldg.
8-11, Kita-Shinagawa 1-chome
Shinagawa-ku, Tokyo, 140-0001 Japan

Phone +81.3.4331.4300
Fax +81.3.4331.4301

Malaysia

Sartorius Malaysia Sdn. Bhd
Lot L3-E-3B, Enterprise 4
Technology Park Malaysia
Bukit Jalil
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622
Fax +60.3.8996.0755

Singapore

Sartorius Singapore Pte. Ltd
1 Science Park Road,
The Capricorn, #05-08A,
Singapore Science Park II
Singapore 117528

Phone +65.6872.3966
Fax +65.6778.2494

South Korea

Sartorius Korea Ltd.
8th Floor, Solid Space B/D,
PanGyoYeok-Ro 220, BunDang-Gu
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700
Fax +82.31.622.5799

Thailand

Sartorius (Thailand) Co. Ltd.
129 Rama 9 Road,
Huaykwang
Bangkok 10310

Phone +66.2643.8361-6
Fax +66.2643.8367



► www.sartorius.com