

PC RECOMMENDED REQUIREMENTS

The PC hosting FLASH & GO must comply with some minimum requirements to support both the Flash software, and the Image Capture hardware and software:

- Pentium IV, 1.7 GHz Processor or higher .
- Windows XP updated with last versions of DirectX and Dot Net utilities or Windows Vista (x32 Edition).
- 1024 Mbytes ram memory or higher, available for the application.
- 500 Mbytes free, after all the operational system updates are installed.
- CD-ROM disk drive
- The unit need:
 - 1 USB 2.0 and 1 RS232
 - 2 USB 2.0

TECHNICAL SPECIFICATIONS

- Reflected Lighting: Ring of 54 white LED.
 - Transmitted Lighting: Two High Flux LED.
 - Background field: Dark or Bright.
 - Imaging device: CMOS with fixed focus lens.
 - Image field: 97x97 mm.
 - Minimum colony size: 0.06 mm.
 - Image capture: True color, 24 bits per pixel.
 - Image resolution: 1536x1536 pixels.
 - Mains: 100/240 V AC, 50/60 Hz
 - Power Consumption: 40W.
 - Operating Temperature: 5°C to 40°C.
 - Operating Relative Humidity: 10% to 90%.
 - Dimensions (WxHxD): 18x36x22 cm.
 - Weight: 5.4 Kg.
- Specifications subject to change without notice

ORDERING INFORMATION

	Cat. No.
FLASH & GO, Automatic colony counter for pour plates and any type of spiral spreading. Complete instrument with centering holder for different Petri dishes size. Software to count colonies in any media. With USB 2.0 high resolution camera to be connected to an external PC. The PC is not included. To run the instrument also monitor, keyboard, and mouse are needed. A notebook with 2 USB 2.0 port can also be used.	6010
Upgrade advanced automatic colony counter	6011
Upgrade advanced inhibition zone measuring	6012
Upgrade user permission administration	6013
Upgrade process selection automation	6014
Plate Handler	4300
Hand for Flash & Go	4888
Base for Flash & Go	6082



Contact us in internet.

To see how the instruments works in a multimedia demonstration, to buy instruments, accessories or consumables on-line visit our web page:

www.iul-inst.com

The IUL instrument owners who are registered in our web will have access to extra services like example of microbiology applications with our instruments, software updates, direct technical service and more.




6146R03 - Mar09

Information

FLASH & GO



www.iul-inst.com

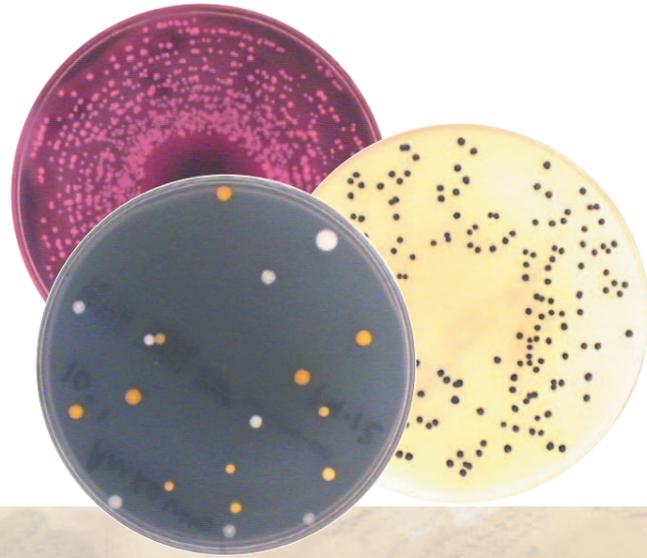


All kinds
of colonies.....

doesn't matter what kind
of colonies are you
counting, the instrument
comes with a high
definition colour camera
which detects colonies of
1/450 inch approx.

.....In different media types can be
counted

The Flash unit has a new illumination
device. The illumination of the plate is
done through a reflection ring. The
background can be black or illuminated
in order to count colonies by transparency.

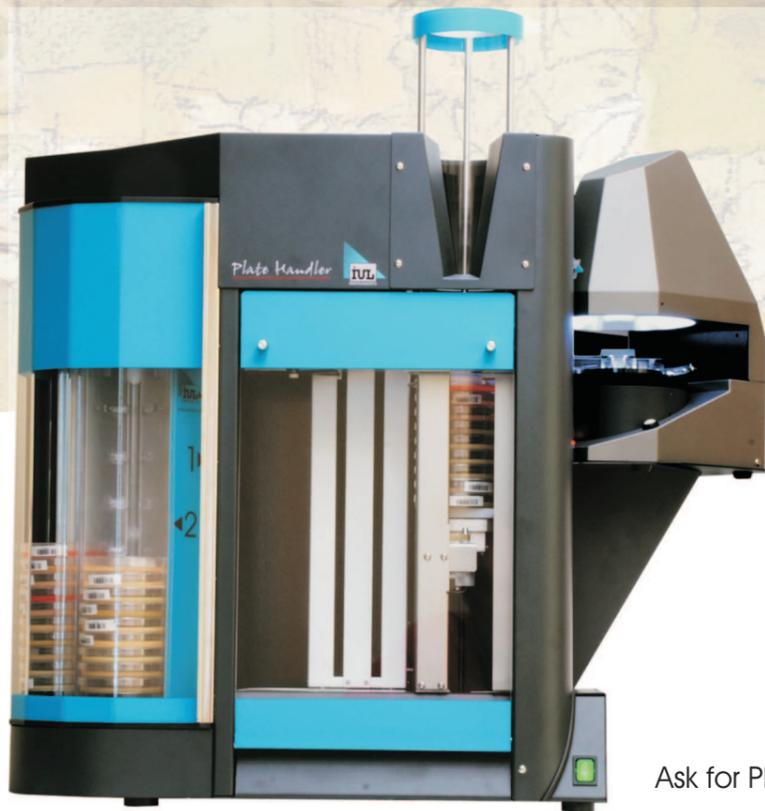

Automation, Connectivity
and use of the
information

Flash & Go is an
instrument to be used in
conjunction with an auto
sampler: the Plate
Handler. All the
information generated by
the Flash such as
recount results or Petri dish
images can be exported
to any software like data
base, spread sheets or
word processors.


Ask for Plate Handler specific leaflet.

The basic software
allows to count up
to six different
types of colonies /
media.

The counting the
analysis of spiral
spreading petri dishes
can be archived by
the basic or upgrade
software.

The upgrade
automated software
is used to define
programs taking in
account colours of
colonies.

A specific software
to measure the size
of inhibitions zones
can be acquired
separately.

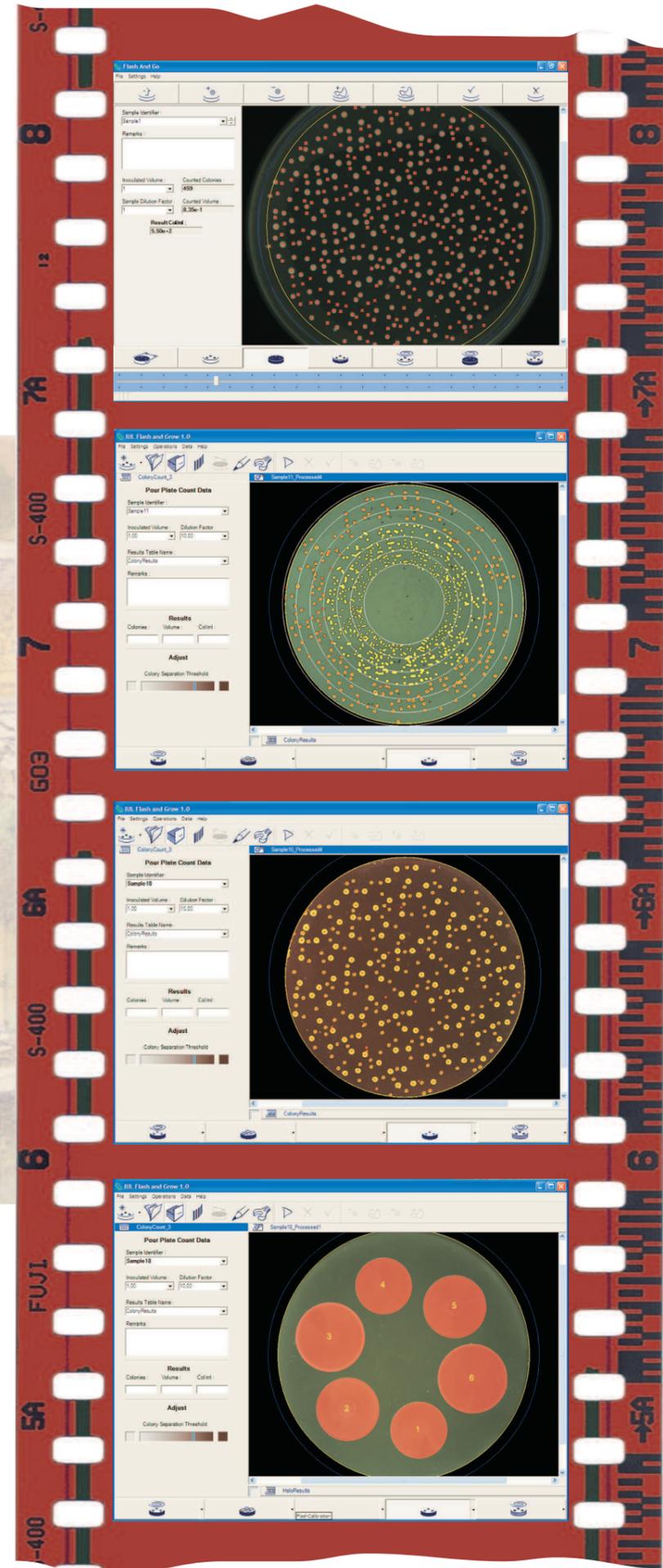

Features and operation